

Web: www.straval.com

Model BPS09i-TC

In-Line Sanitary Tri-Clamp Flange Back Pressure/Bypass Valve



-In-Line, Self-Draining Design

- -Diaphragm or Piston and Spring Operated
- -1/2"-3" Sanitary Tri-Clamp Connections
- -Inlet Pressures Up To 150 PSIG (10.3 BAR), Set Pressure Range From 0-80 PSIG (0-6.9 BAR)
- -In Line Serviceable

-Internal 20 RA Surface Finish Standard Constructed of 316L Stainless Steel

-ANSI Class VI Seat Leakage

Features

- Pressure-Containing Parts made from solid bar stock materials unlike castings which have wall thickness variations.
- Body and Trim: Standard material is type 316L stainless steel mechanically polished to a uniform 20 RA surface finish. Electropolishing for enhanced surface finish is offered at an additional cost.
- Diaphragm or Piston Sensing Mechanism: Diaphragm models offered only in ½"-2" sizes. Diaphragm sensed back pressure valves offer superior sensitivity to increases in process line pressures as opposed to piston sensed valves. However, piston sensed designs are capable of much larger flow rates which are offered in 2-1/2" and 3" connections.
- Teflon-Viton Composite Reinforced Diaphragm: Viton backed diaphragm with a Teflon film interacting on the wetted side offers adequate corrosion resistance to a variety of fluids and gases including steam with an allowable temperature of 350°F (176°C). Diaphragms with Buna backing are available for lower temperature applications. Flexible all Buna or all Viton diaphragms are available as well and are recommended for low pressure applications.
- Spring Chamber and Adjusting Spring: These parts are non-wetted, not in contact with the fluid or gas, and standard in 316L stainless steel.
- In-Line Porting: In line Sanitary Tri-Clamp connections are standard. Mounting in the horizontal upright position is highly recommended as it is the only position that drains water away from the valve trim, otherwise the valve can operate in any orientation.
- Inline Serviceable: Capable of being serviced inline. The process lines must be depressurized before any servicing of this product. All trim components can be replaced, inspected, or cleaned by removing the sanitary clamp.
- Self-Draining: The inclined toroidal geometry connecting the inlet to the flow chamber allows the valve to self-drain.
- 300 Series Stainless Steel V-Band Clamp: Allows for quick assembly and disassembly. Clamp is rated for 800 psi.

Applications

This product is used for sanitary backpressure regulation. A few typical applications include maintaining constant pump discharge pressures, bypassing excessive pressures from various types of process equipment, and any application where a constant pressure must be maintained in a sanitary process or piping system. This product is designed for clean, filtered or strained, non-abrasive, liquids and gases including but not limited to water, steam, non-fluorinated chemicals, inert gasses and fluids compatible with the selected elastomers and valve materials. A strainer or filter with the appropriate material and pressure rating can be purchased from STRAVAL. High temperature applications can be satisfied with careful selection of materials (contact factory).

Options

- ◆ Connection Sizes: ½"-3" tri-clamp sanitary ferrules typical. See BPS09i-FLG or BPS09i-NPT for industrial connections.
- Diaphragm Sensing (½"-2" Only)
 Viton Backing with PTFE Film (wetted side), Buna Backing with PTFE Film (wetted side), Only Buna, Only Viton
 - Seat: PTFE, Buna, Viton, Metal
- ◆ Piston Sensing (2-½"-3" Only)
 - Body and Spring Chamber Seal: EPDM, Buna, PTFE, Viton, Kalrez
 - Piston Seal: EPDM, Buna, PTFE, Viton, Kalrez
 - Seat: EPDM, Buna, PTFE, Viton, Kalrez
- Surface Finish: 20 RA mechanical polish, Electropolished

*Price Varies with Materials of Construction, Size, and Selected Options/Alternatives


Principle of Operation

This is a direct acting valve with an adjustable spring operating against a diaphragm or piston subjected to the inlet

pressure of the valve. Increasing the spring compression by turning the adjusting screw clockwise, increases the system or line pressure to be maintained. Reducing the spring compression by turning the adjusting screw counter-clockwise will reduce the system or line pressure to be maintained. An increase in line pressure beyond the set point (inlet pressure that exceeds the compressive force of the spring) will cause the seat or diaphragm to open permitting flow to the outlet. This bypasses any excess line pressure through the outlet of the valve

Valve Type	Straval Identification Number	Porting	Connection Size	Connection Type	Sensing Mechanism
BPS-Back Pressure Regulator	09	i-In-Line	05: %" 07: %" 10: 1" 15: 1-%" 20: 2" 25: 2-%" 30: 3"	San. Tri-Clamp: TC	D-Diaphragm D-Diaphragm D-Diaphragm D-Diaphragm D-Diaphragm P-Piston P-Piston
	E	3PS 09	i-10TC	P	PPISON

DIAPHRAGM SENSED

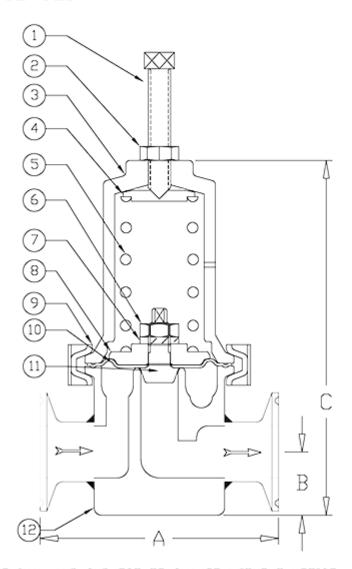

ALL DIMENSIONS NOMINAL AND SUBJECT TO CHANGE WITHOUT NOTICE CONTACT FACTORY FOR CERTIFIED DIMENSIONAL DRAWINGS **DIMENSION C IS DEPENDANT ON SPRING RANGE, CONTACT PLANT

MATERIAL LIST			DIMENSIONS (inches)				
& SPECIFICATION			SIZE	cv	А	В	С
1	PART	MATERIAL COLUMN SECONOMICS	1/2	1.85	4'	13/16*	**
	ADJUSTING SCREW	304, 316L STAINLESS STEEL	17.6	1.00		107.10	
S	JAM NUT	304, 316L STAINLESS STEEL	3/4	3.1	4'	15/16*	××
3	SPRING CHAMBER	303, 316L STAINLESS STEEL	1	10.3	6'	1"	××
4	SPRING PUSHER	303, 316L STAINLESS STEEL	1-1/2	12.2	6'	1 1/4"	**
5	SPRING	304, 316L STAINLESS STEEL	s	12.2	6*	1 1/2*	××
6	NUT	304, 316L STAINLESS STEEL					
7	LOCK WASHER	304, 316L STAINLESS STEEL	PH 973-345- FX 973-340-		-337	13/8 VM	NES
8	SPRING FOLLOWER	303, 316L STAINLESS STEEL	ELAVOCO PAR	WE	STRAI	NERS VA	77,
9	V-BAND CLAMP	304 STAINLESS STEEL	NATL TARRESTOR	rine Par	ela Danceria	re/Bypass	Value
10	DIAPHRAGM	TFE/Viton,TFE/Buna,Buna,Vitor		DN-U	LINE SANITARY		44146
11	DIAPHRAGM SCREW	303, 316L STAINLESS STEEL	BATE 08/15/2008	29/5 AO		TO D	REV.
12	BODY	303, 316L STAINLESS STEEL	DRVW AS] #	3PS09i-	- I C - D	Α.


Material List and Specification

Diaphragm Sensed

#	Part	Materials		
1.	Adjusting Screw	304, 316L Stainless Steel		
2.	Jam Nut	304, 316L Stainless Steel		
3.	Spring Chamber	303, 316L Stainless Steel		
4.	Spring Pusher	303, 316L Stainless Steel		
5.	Spring	304, 316L Stainless Steel		
6.	Nut	304, 316L Stainless Steel		
7.	Lock Washer	304, 316L Stainless Steel		
8.	Spring Follower	303, 316L Stainless Steel		
9.	V-Band Clamp	304 Stainless Steel		
10.	Diaphragm	TFE/Vitorn, TFE/Buna, Buna, Vito		
11.	Diaphragm Screw	304, 316L Stainless Steel		
12.	Body	303, 316L Stainless Steel		

Dimensions (inch)						
Size	cv	A	В	C		
1/2	1.85	4"	13/16"	**		
3/4	3.1	4"	15/16"	**		
1	10.3	6"	1"	**		
1-1/2	12.2	6"	1 1/4"	**		
2	12.2	6"	1 1/2"	**		

Note: Dimensions are approximate and are subject to change without notice. Request certified dimensions before final product installation.

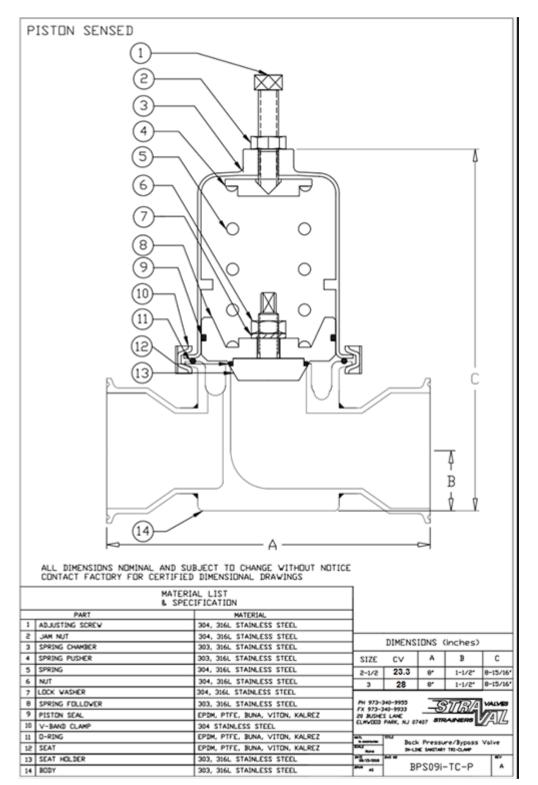

Material List and Specification

Piston Sensed

#	Part	Materials		
1.	Adjusting Screw	304, 316L Stainless Steel		
2.	Jam Nut	304, 316L Stainless Steel		
3.	Spring Chamber	303, 316L Stainless Steel		
		303, 316L Stainless Steel		
5.	Spring	304, 316L Stainless Steel		
6.	Nut	304, 316L Stainless Steel		

7.	Lock Washer	304, 316L Stainless Steel		
8.	Spring Follower	303, 316L Stainless Steel		
9.	Piston Seal	EPDM, PTFE, Buna, Viton, Kalrez		
10.	V-Band Clamp	304 Stainless Steel		
11.	O-Ring	EPDM, PTFE, Buna, Viton, Kalrez		
12.	Seat	EPDM, PTFE, Buna, Viton, Kalrez		
13.	Seat Holder	303, 316L Stainless Steel		
14.	Body	303, 316L Stainless Steel		

Dimensions (inch)					
Size	cv	A	В	C	
2-1/2	23.3	8"	1-1/2"	8-15/16"	
3	28	8"	1-1/2"	8-15/16"	

Note: Dimensions are approximate and are subject to change without notice. Request certified dimensions before final product installation.

1/2" BPS09i-05TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from:2-80 psig (0.138-5.52 barg) Select spring from pricing page

3/4" BPS09i-07TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from: 2-80 psig (0.138-5.52 barg) Select spring from pricing page

1" BPS09i-10TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from: 2-80 psig (0.138-5.52 barg) Select spring from pricing page

1 1/2" BPS09i-15TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from: 2-80 psig (0.138-5.52 barg) Select spring from pricing page

2" BPS09i-20TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from:2-80 psig (0.138-5.52 barg) Select spring from pricing page

2 1/2" BPS09i-25TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from: 2-80 psig (0.138-5.52 barg) Select spring from pricing page

3" BPS09i-30TC

Sanitary Tri-Clamp Connections 150 psi (~10 bar) rating

Multiple Spring Ranges from: 2-80 psig (0.138-5.52 barg) Select spring from pricing page

The spring ranges listed above are not achievable with one spring, but are compressed to show overall product capability. Select a specific spring range in the pricing pages or specify a set pressure when ordering.